

GeneSpin Srl Via Friuli, 51 -20135 Milano P.IVA e C.F. 04520270960 administration@pec.genespin.com info@genespin.com

## > Blasticidin S HCI

cat. no.	amount	note
STS-BLA25	2 <b>5</b> mg	powder
STS-BLA21	100mg	powder

Blasticidin S is a peptidyl nucleoside antibiotic isolated from the culture broth of Streptomyces griseochromogenes. Blasticidin S specifically inhibits protein synthesis in both prokaryotes and eukaryotes through inhibition of peptide-bond formation in the ribosomal machinery.

Blasticidin S allows the selection and maintenance of cells expressing the blasticidin-resistance gene. Three blasticidin-resistance genes have been cloned and sequenced: an acetyl transferase gene, bls, from Streptoverticillum sp. JCM 4673, a blasticidin producer strain, and two deaminase genes, bsr, from Bacillus cereus and BSD from Aspergillus terreus. Both genes, bsr and BSD, are used as dominant selectable markers for transformation experiments in mammalian and plant cells. Although Blasticidin S was developed as a selection agent for mammalian cells, the bsr and BSD resistance genes can also be used in E. coli.

Blasticidin S resistant transformants can be selected in low salt LB agar medium, pH 8, supplemented with 100  $\mu$ g/ml blasticidin S. The working concentrations of blasticidin S for mammalian cell lines varie from 3 to 50  $\mu$ g/ml. In a starting experiment we recommend to determine optimal concentrations of antibiotic required to kill your host cell line. After treatment, cell death occurs rapidly allowing the selection of transfected cells with plasmids carrying the BSD or bsr genes in as little as 7 days post-transfection.

Formulation  $C_{17}H_{26}N_8O_5$ Molecular Weight 458.9

CAS No. 3513-03-9 Purity: >98.0%

Solubility Solubility in water..

Storage Store dry at 0-5°C.

FOR RESEARCH USE ONLY



## > Blasticidin S HCI Solution

antibiotic

cat. no.	amount	note
STS-BLA20	20mg	solution
STS-BLA21	100mg	solution

Blasticidin S is a peptidyl nucleoside antibiotic isolated from the culture broth of Streptomyces griseochromogenes. Blasticidin S specifically inhibits protein synthesis in both prokaryotes and eukaryotes through inhibition of peptide-bond formation in the ribosomal machinery.

• DATA SHEET

Blasticidin S allows the selection and maintenance of cells expressing the blasticidin-resistance gene. Three blasticidin-resistance genes have been cloned and sequenced: an acetyl transferase gene, bls, from Streptoverticillum sp. JCM 4673, a blasticidin producer strain, and two deaminase genes, bsr, from Bacillus cereus and BSD from Aspergillus terreus. Both genes, bsr and BSD, are used as dominant selectable markers for transformation experiments in mammalian and plant cells. Although Blasticidin S was developed as a selection agent for mammalian cells, the bsr and BSD resistance genes can also be used in E. coli.

Blasticidin S resistant transformants can be selected in low salt LB agar medium, pH 8, supplemented with 100  $\mu$ g/ml blasticidin S. The working concentrations of blasticidin S for mammalian cell lines varie from 3 to 50  $\mu$ g/ml. In a starting experiment we recommend to determine optimal concentrations of antibiotic required to kill your host cell line. After treatment, cell death occurs rapidly allowing the selection of transfected cells with plasmids carrying the BSD or bsr genes in as little as 7 days post-transfection.

Formulation C<sub>17</sub>H<sub>26</sub>N<sub>8</sub>O<sub>5</sub> Molecular Weight 458,9

Solubility Supplied as a 10mg/ml sterile solution in HEPES buffer. Soluble in water and slightly alkaline or acid solution.

Storage Store dry at -20°C.