



DATA SHEET

info@genespin.com - www.genespin.com

GeneSpin Srl Via Friuli, 51 -20135 Milano
P.IVA e C.F. 04520270960
administration@pec.genespin.com
info@genespin.com

> K Large DNA Ladder

A unique combination of a number of proprietary plasmids digested with appropriate restriction enzymes and PCR products to yield 19 fragments, suitable for use as molecular weight standards for agarose gel electrophoresis. The DNA includes fragments ranging from 100-10,000 base pairs. The 500, 1.5K and 3K bands have increased intensity to serve as reference points. The approximate mass of DNA in each band is provided (0.86 µg a load) for approximating the mass of DNA in comparably intense samples of similar size.

cat. no.	amount	note	DNA fragments
STS-KL	500ul	range: 100bp to 10Kb	0.1, 0.2, 0.3, 0.4, 0.5, 0.6, 0.7, 0.8, 0.9, 1.0, 1.5, 2.5, 3.0, 4.0, 5.0, 6.0, 8.0, 10kb

SOURCE:

PCR products and double-stranded DNA digested with appropriate restriction enzymes, are phenol extracted and equilibrated to 10mM Tris-HCl (pH 8.0) and 1mM EDTA.

RANGE and BANDS

100-10000bp
19 bands

RECOMMENDED LOAD: 5ul/well


Containing orange G & xylene cyanol FF as tracking dyes.

STORE:

Store at 4°C for 12 month.

Under appropriate Storage Conditions, quality is guaranteed for 24 months at -20°C .

StoS KL Ladder

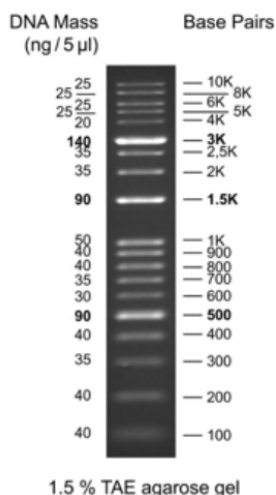

FOR RESEARCH USE ONLY